

Chem CD

The Formaldehyde Replacement



Chem – CD is a unique product specifically aimed at replacing Formaldehyde and other methods for sterilisation of BSC's and Isolators.

Easy to use, easy to store, cost effective and most importantly a highly effective way of ensuring a sterile environment.

*Chem CD comes in 3 parts, A, B and N, all of which are pre-measured sachets for ease of use. The gas produced leaves **no residues** and is **non-corrosive** to protect the inside of the target chamber.*

For more information contact:

enquiries@primattec.co.uk

or call 01722 33 66 33



TO USE:

1. **Simply mix parts A+B with water.**
2. **Seal your BSC / Isolator.**
3. **Leave to produce gas and sit for 85 minutes.**
4. **Start carbon scrubber system**
5. **When chamber is empty mix Part N with water, open the chamber and pour into parts A+B.**
6. **You now have a Neutral salt solution to pour down the**